


## Body Belt Incubator – Users' Guide

The body belt incubator is an easy to use, cost effective and a reliable technique to be used in field for incubating bacterial plates. Each body belt incubator has at least 8 pockets to carry bacterial plates, and a pocket for thermometer.

- 1) After the test, close the lid of the plates correctly and place the plates in the body belt incubator pockets.
- 2) Place the thermometer in its respective pocket.
- 3) Wear the belt around the belly for 24 to 48 hours.

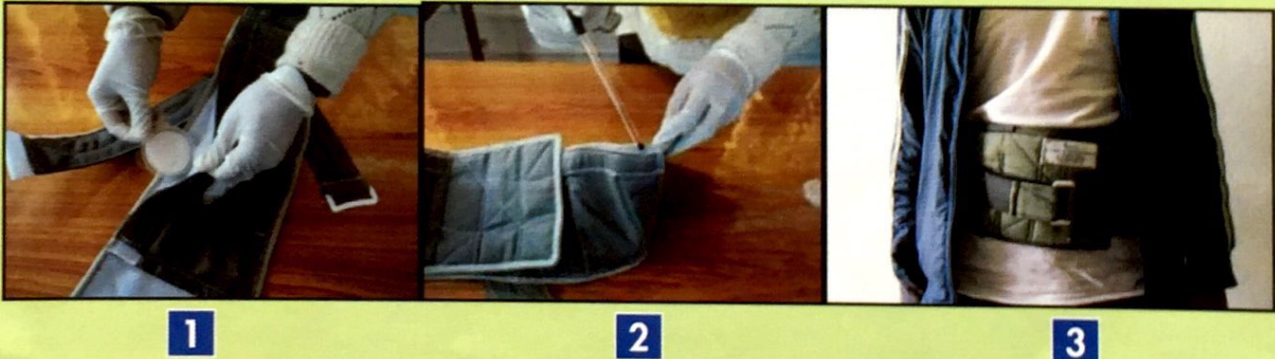

### Things to note for better performance of the Body Belt Incubator:

- Body Belt Incubator **should be** worn regularly for 24 hours.
- Body Belt Incubator **should not be** worn over thick clothes. It is advisable to have the belt in body contact as far as possible; wear it over thin clothes.
- The temperature **must be** regularly monitored using the thermometer.





संस्थापक र जनसंचालक संस्था (एनपीओ)  
 पते: बन्स नं: ४१०२, बानेश्वर मार्ग, बापामार्ग, नयाँ बानेश्वर, काठमाडौं ।  
 फोन नं: ०१-४२४४४४४, ४२४४४४४, फ्याक्स: ०१-४२४४४४४  
 ईमेल: enpho@enpho.org / वेब साइट: www.enpho.org



Eco Concern Pvt. Ltd.  
 New Baneshwor, P. O. Box: 3456  
 Kathmandu, Nepal  
 Phone: +977-1-5244301

## बडि बेल्ट इन्क्यूबेटर : प्रयोगकर्ता निर्देशिका

बडि बेल्ट इन्क्यूबेटर फिल्ट्रमा ब्याक्टेरीया इन्क्यूबेशन गर्न एक सजिलो, लागत प्रभावकारी र विश्वसनीय प्रविधि हो । प्रत्येक बडि बेल्ट इन्क्यूबेटरमा जिवाणु प्लेट राख्न कम्तिमा ८ वटा पकेट र थर्मोमिटर राख्न १ पकेट हुन्छ ।

- १) टेष्ट गरिसकेपछि, राम्रोसँग जीवाणु प्लेट बन्द गर्ने र बडि बेल्ट इन्क्यूबेटर पकेटमा जिवाणु प्लेट राख्ने ।
- २) थर्मोमिटर निर्धारित पकेटमा राख्ने ।
- ३) बडि बेल्ट इन्क्यूबेटर कम्मरमा २४ देखि ४८ घण्टा सम्म लगाउने ।



१

२

३

बडि बेल्ट इन्क्यूबेटरको प्रभावकारीता को लागि ध्यान दिनुपर्ने कुराहरू :

- बडि बेल्ट इन्क्यूबेटर कम्मरमा लगातर २४ घण्टा लगाउनु पर्छ ।
- बडि बेल्ट इन्क्यूबेटर बाक्लो कपडामाथि लगाउनु हुँदैन ।
- बडि बेल्ट इन्क्यूबेटर शरिरसँग सिधा सम्पर्क हुने गरि वा पातलो कपडामाथि लगाउनु पर्छ ।
- थर्मोमिटरको प्रयोग गरी तापक्रमको लगातर अनुगमन गर्नुपर्छ ।